

# User's Guide



*Shop online at*

**omega.com®**

Ω OMEGA®

*omega.com*

*e-mail: info@omega.com*

*For latest product manuals:*

*omegamanual.info*

**ISO 9001**  
CERTIFIED  
CORPORATE QUALITY

STAMFORD, CT

**ISO 9001**  
CERTIFIED  
CORPORATE QUALITY

MANCHESTER, UK

## TCY 20, 25, 30 & 48 Thermal Cyclers



**OMEGAnet® Online Service**  
**omega.com**

**Internet e-mail**  
**info@omega.com**

### **Servicing North America:**

**U.S.A.:**  
ISO 9001 Certified

Omega Engineering, Inc., One Omega Drive, P.O. Box 4047  
Stamford, CT 06907-0047 USA  
Toll Free: 1-800-826-6342  
FAX: (203) 359-7700

TEL: (203) 359-1660  
e-mail: info@omega.com

**Canada:**

976 Bergar  
Laval (Quebec), Canada H7L 5A1  
Toll-Free: 1-800-826-6342  
FAX: (514) 856-6886

TEL: (514) 856-6928  
e-mail: info@omega.ca

### **For immediate technical or application assistance:**

**U.S.A. and Canada:** Sales Service: 1-800-826-6342/1-800-TC-OMEGA®  
Customer Service: 1-800-622-2378/1-800-622-BEST®  
Engineering Service: 1-800-872-9436/1-800-USA-WHEN®

**Mexico:**

En Español: 001 (203) 359-7803  
info@omega.com.mx

FAX: (001) 203-359-7807  
e-mail: espanol@omega.com

### **Servicing Europe:**

**Benelux:**

Managed by the United Kingdom Office  
Toll-Free: 0800 099 3344  
FAX: +31 20 643 46 43

TEL: +31 20 347 21 21  
e-mail: sales@omega.nl

**Czech Republic:**

Frystatska 184  
733 01 Karviná, Czech Republic  
Toll-Free: 0800-1-66342  
FAX: +420-59-6311114

TEL: +420-59-6311899  
e-mail: info@omegashop.cz

**France:**

Managed by the United Kingdom Office  
Toll-Free: 0800 466 342  
FAX: +33 (0) 130 57 54 27

TEL: +33 (0) 161 37 29 00  
e-mail: sales@omega.fr

**Germany/Austria:**

Daimlerstrasse 26  
D-75392 Deckenpfronn, Germany  
Toll-Free: 0 800 6397678  
FAX: +49 (0) 7056 9398-29

TEL: +49 (0) 7059 9398-0  
e-mail: info@omega.de

**United Kingdom:**  
ISO 9001 Certified

OMEGA Engineering Ltd.  
One Omega Drive, River Bend Technology Centre, Northbank  
Irlam, Manchester M44 5BD England  
Toll-Free: 0800-488-488  
FAX: +44 (0)161 777-6622

TEL: +44 (0)161 777-6611  
e-mail: sales@omega.co.uk

It is the policy of OMEGA Engineering, Inc. to comply with all worldwide safety and EMC/EMI regulations that apply. OMEGA is constantly pursuing certification of its products to the European New Approach Directives. OMEGA will add the CE mark to every appropriate device upon certification.

The information contained in this document is believed to be correct, but OMEGA accepts no liability for any errors it contains, and reserves the right to alter specifications without notice.

**WARNING:** These products are not designed for use in, and should not be used for, human applications.

## Contents

<b>1. Introduction .....</b>	<b>4</b>
<b>2. Safety and installation.....</b>	<b>4</b>
<b>3. Cyclor TCY specification.....</b>	<b>7</b>
<b>4. User operation and consumables .....</b>	<b>8</b>
Heated lid .....	8
Tubes or reaction vessels.....	8
0.5ml Tubes.....	9
Switching on .....	10
<b>5. Front panel controls .....</b>	<b>11</b>
The LCD Display .....	11
Indicators.....	11
The keys.....	12
<b>6. Programming the Cyclor TCY.....</b>	<b>13</b>
Programming Options.....	15
Option 1: Creating programs using templates.....	19
Option 2: Creating a new program.....	20
Example programs .....	20
Programming.....	20
To program increment or decrement temperature or time. ....	24
To run a program.....	27
<b>7. Information .....</b>	<b>29</b>
<b>8. Configuration .....</b>	<b>29</b>
<b>9. Notes on transferring protocols from other thermal cyclers.....</b>	<b>32</b>
<b>10. User maintenance .....</b>	<b>33</b>
Cleaning your Cyclor TCY .....	33
Heater Block removal .....	33
<b>11. Some frequently asked questions .....</b>	<b>34</b>
<b>12. Additional information .....</b>	<b>35</b>
Fault Finding.....	35
Fuses .....	35
The heated lid over-temperature cut-out.....	35
Insulation testing .....	35
Interchangeable blocks.....	35

## 1. Introduction

The Cyclor TCY provides the researcher with the means of accurately controlling the temperature profile of samples. It has many scientific applications, including DNA amplification and sequencing. The Cyclor TCY can cycle samples between 4°C and 99°C.

The Cyclor TCY is programmed by means of an integral keypad and LCD display. A program, which can be recalled from memory, consists of:

- a series of specified temperatures in °C
- the times for which each specified temperature will be held (Hold Times)
- the desired heating or cooling rates, in °C/sec, between each specified temperature
- optional instructions as to whether the times and/or the temperatures are to increase or decrease when a cycle is repeated.

The memory can store up to 80 programs.

Cyclor TCY is the most economic personal thermal cyclor. There are four types of blocks can be chosen for this model: 0.2ml X 25 and 0.5ml X 20 wells blocks which have less sample compartment but with higher ramping rate; 0.2 ml X 48 and 0.5ml X 60 wells blocks which have more wells but less ramping rate.

Before using the TCY units, make sure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment contact Omega Flow Engineering for assistance.

## 2. Safety and installation

### **Operator Safety**

Please read this manual carefully before using the thermal cyclor. If the equipment is not used in the manner described in this manual the protection provided by the equipment might be impaired.

All Omega Engineering, Inc. instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out. If a safety problem should be encountered then switch off the unit at the power source and remove the plug from the electricity supply.

### **Warning**

HIGH TEMPERATURES ARE DANGEROUS as they can cause serious burns to operators and ignite combustible material. Users should be aware of the following potential hazards:



- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS
- DO NOT use combustible substances near hot objects
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases
- DO NOT place any liquid directly into the instrument, always use the appropriate plastic consumables

After heating samples, remember that parts of the instrument, namely the lid and block and any associated accessories may be very hot.

## ***Unpacking***

---

When unpacking please ensure that the following have been removed from the packaging:

- Thermal cyclers
- Power cable

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Omega Engineering, Inc. accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

Please note that the screen of the thermal cycler is easily damaged by sharp objects such as pens, pencils and fingernails. This type of damage will be considered as misuse and invalidate the guarantee for this component.

## ***Installation***

---

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the power electricity supply.

1. All Omega Engineering, Inc. instruments are supplied with a power cable; this may be integral or plug-in.
2. Before connecting the instrument to the power electricity supply, check the voltage against the rating plate (located on the back of the unit). Ensure that the voltage selector switch (located above the On/Off switch) is set to the appropriate voltage for the local supply. Please note that the unit must be grounded to ensure proper electrical safety. Connect the power cable to a suitable plug according to the table below.

Connection	220/240V, 50/60Hz Supply	110V/120V Supply, 50/60Hz
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

3. Units showing 230V, 50/60Hz on the rating plate also operate between 210 and 260V; units with 120V, 50/60Hz also operate between 100 and 130V and units marked with 100V will operate between 90 and 110V. In all cases the heating rate will degrade by approximately 8% at the extremes of the voltage range.
4. Plug the power cable into the socket on the back of the instrument.
5. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction. Position the instrument with a minimum distance all round of 200mm from walls or other items and between each unit if multiple units are being used to permit efficient air flow for each instrument.
6. Switch on the instrument:

I      Power Switch On

O      Power Switch Off

### **Replacement cable**

Should the power lead need replacement, a cable of 1mm<sup>2</sup> of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

### ***Working Conditions***

---

The thermal cycler is designed operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Relative humidity not exceeding 95%
- Power supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of 10°C to 30°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

### **3. Cyclor TCY Specification**

#### ***Temperature***

---

Temperature range	4°C to 99°C
Temperature set point precision	0.1°C
Block uniformity (over full range)	± 0.5°C
Temperature accuracy	± 0.5°C
End of program cooling below ambient	Yes

#### ***Heating/Cooling Rate***

---

Maximum heating rate	4.5°C/s
Minimum heating rate	0.1°C/s
Maximum cooling rate	3.0°C/s
Minimum cooling rate	0.1°C/s

#### ***Heated Lid***

---

Heated lid enable/disable	Yes
Selectable heated lid temperature	100°C to 115°C
Over-temperature cut-out	Fixed at 145°C
Temperature sensor	Thermistor
Heater type	Etched foil
Heater power	33W
Warm up time	4 min approximately (ambient to 105°C)

The heated lid will only come on if the set temperature is above 35°C

#### ***Programming***

---

Program naming using alpha keys	Yes
Program password protection	Yes
Number of programs (3 step)	37
Maximum number of steps per stage	35
Maximum number of stages per program	25
Minimum hold time	1 sec
Maximum hold time	18 hrs
Incremental/decremental temperature	Yes
Incremental/decremental step hold time	Yes
Programmable ramp rate (heat or cool)	Variable in 0.1°C/sec steps
Programmable ramp rate resolution	0.1°C

#### ***Running Programs***

---

Pause button	Yes
Stop button	Yes
End of program alarm	Yes
Auto restart on power fail	Yes

#### ***Interconnectivity***

---

Connection to PC control program	Yes
Software updates from web-site	Yes

### **Dimensions**

Height	185mm
Width	185mm
Length	330mm

### **Power**

Power consumption	230W
-------------------	------

Units marked 100V will work at any voltage between 90 and 110V.  
The performance will vary and will not necessarily meet the above typical specification at the extremes of voltage.

## **4. User operation and consumables**

### **Heated lid**

To release the heated lid, lift the lid catch and the lid will open. To close the lid, push down gently until the catch clicks in place. **Do not close the lid without tubes being fitted in the block.**



### **Tubes or reaction vessels**

Omega Engineering, Inc. does not recommend any specific tube or reaction vessel other than those described in this manual. We recommend using reaction volumes between 20 and 200 $\mu$ l. The tubes must withstand a pressure of 1 atmosphere at 100°C. Any vessel must be able to withstand the temperatures you are using without any danger of them deforming to the point where they fracture.

To test your tubes, put 25 $\mu$ l of water into each of 5 tubes and subject them to a typical thermal cycling protocol. At the end of the program, measure the volume remaining using a micropipette. A loss of more than 1 or 2 $\mu$ l indicates a vapour leak.




The amount of volume loss you observe and the change in reactant concentrations you can tolerate determine the minimum volume that can be used. Typical volume losses of 1µl in 30 cycles allow the use of samples of 20µl or less.

During the final cool-down, a ring of condensation may form above the liquid level but below the top of the sample block. This is normal and does not effect the reaction.

### **0.5ml Tubes**


The heated lid mechanism of the Cycler TCY with the 0.5ml block is optimised for flat top tubes. In order to fit the maximum number of tubes in the block, they need to be placed so that all the hinges are in the same relative orientation. An example is illustrated below.

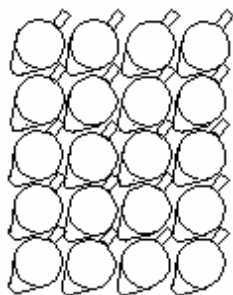


**Place the tubes in the block in the same orientation in order to achieve the maximum capacity.**

**Always make sure the pressure on the heated lid is balanced by spreading the tubes across the block or inserting some dummy tubes if necessary.**

## Switching on

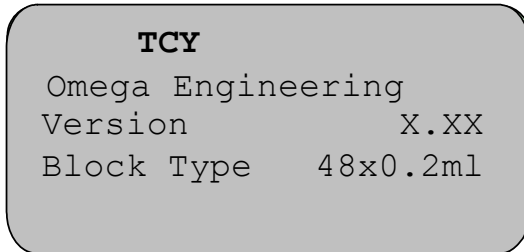

When you switch on the unit a screen similar to the one shown here will appear briefly. This indicates the software version and block type installed in the unit.

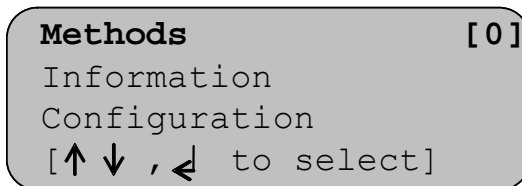


When the initial screen has disappeared the screen shown here will automatically appear. There are three options:


1. 'Programs' for creating and running programs;
2. 'Information' for obtaining information about the unit e.g. serial number;
3. 'Configuration' for setting up program defaults and settings.

Each of these is explained in detail in section 6.


## 5. Front panel controls

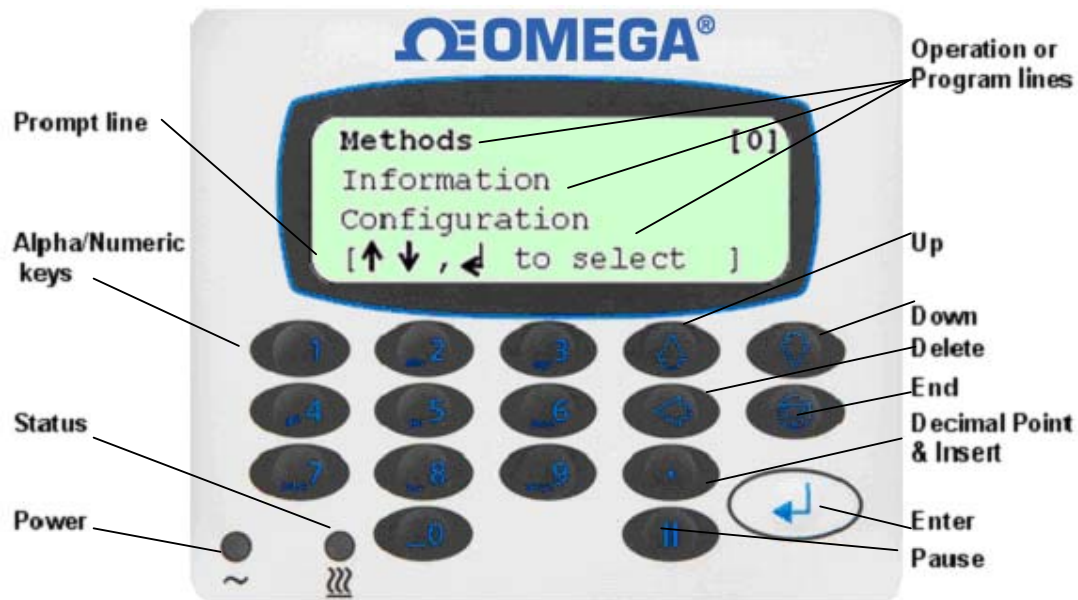

### The LCD Display

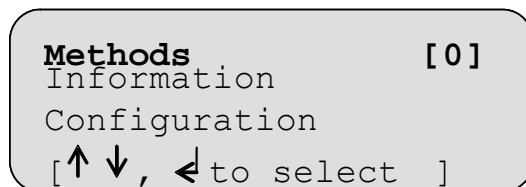

The CYCLER TCY has a four line LCD display. The top three lines are information or program lines, while the bottom line is a prompt line. The active line flashes on screen and is shown in bold in this manual.

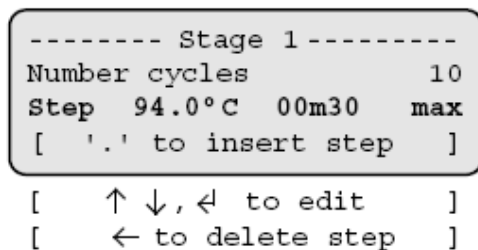

The bottom line may change or flash between two or three different prompts depending upon which program line is active. An example of a prompt line which flashes between three different prompts is shown here.

### Indicators



**Power:** This indicator is lit when there is power to the CYCLER TCY.



**Status:** This indicator is lit when the CYCLER TCY is running a program. It is off when the unit is stopped. It will flash and beep slowly when the unit is paused. It will flash after a five second quick sequence of beeps when the unit reaches its final hold temperature.

## The keys

*The End, Finish or Exit key*



The “End” or “Finish” key is used to finish editing programs or to exit a program. It is also used to exit a sequence of key operations.

*The Up arrow key*



The “Up arrow” key is used for scrolling up through the lines on the screen. It is also used to increase certain fields when editing.

*The Down arrow key*



The “Down arrow” key is used for scrolling down through the lines on the screen. It is also used to decrease certain fields when editing.

*The Decimal Point and Insert key*



The “Decimal Point” key is used when editing numeric fields. It is also used as an Insert key during program editing to insert additional steps and other field sets.

*The Delete key*



The “Delete” key is used during program editing to delete steps and stages.

*The Pause key*



The “Pause” key is used during program editing to insert a pause after a stage or to insert increment and decrement times and temperatures. It is also used to pause a program when it is running.

*The Enter or Accept key*



The “Enter” key is used to enter an operation or the edit mode. It is also used to ‘Accept’ a field or program change.

*The End, Finish or Exit key*



The “End” or “Finish” key is used to finish editing programs or to exit a program. It is also used to exit a sequence of key operations.

*The Up arrow key*



The “Up arrow” key is used for scrolling up through the lines on the screen. It is also used to increase certain fields when editing.

*The Down arrow key*



The “Down arrow” key is used for scrolling down through the lines on the screen. It is also used to decrease certain fields when editing.

*The Decimal Point and Insert key*



The “Decimal Point” key is used when editing numeric fields. It is also used as an Insert key during program editing to insert additional steps and other field sets.

*The Delete key*



The “Delete” key is used during program editing to delete steps and stages.

*The Pause key*



The “Pause” key is used during program editing to insert a pause after a stage or to insert increment and decrement times and temperatures. It is also used to pause a program when it is running.

*The Enter or Accept key*



The “Enter” key is used to enter an operation or the edit mode. It is also used to ‘Accept’ a field or program change.












*The Alpha/Numeric keys*

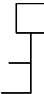

The “Alpha/Numeric” keys are used to enter numbers and letters during programming. If it is a numeric field then numbers will be inserted; if it is an alpha field then letters will be inserted. Each key will cycle through the letters shown on the key (0 will insert a space). To insert a number, press and hold the appropriate key.

## **6. Programming the CYCLER**

The table below illustrates the symbols which appear on the LCD display and the relevant keypad button to which they refer.

	In the prompt line is equivalent to the “End” key.	
	On a program line means the program is locked; it can be copied and then edited but not edited directly.	
	Up arrow	
	Down arrow	
	Enter	
	Pause	

The following terms are used when programming the CYCLER TCY:

<b>Stage</b>	A ‘Stage’ is the name given to the main repeated part of a program containing the ‘Steps’ to be repeated. A program can contain a maximum of 25 stages although typically only one or two are used.
<b>Cycle</b>	A ‘Cycle’ is the number of repetitions of a ‘Stage’
<b>Step</b>	A ‘Step’ is a user defined program line containing a temperature, a hold time and ramp rate. Typically three steps are created within one ‘Stage’.

```

Methods [ 0 ]
Information
Configuration
[ ↑ ↓ , ↵ to select ]

```

To enter the programming section of the Cyclor TCY software, highlight ‘methods’ using the “Up Arrow” or “Down Arrow” keys  
 ↑ ↓  
 Press “Enter” ↵

```

* *   New Methods   * * [ ]
2 STEP               5
3 STEP               5
[ ↑ ↓ , ↵ to select ]

```

This screen lists all the programs saved on the unit. Use the “Up Arrow” and “Down Arrow” keys ↑ ↓ to scroll through the list.

The number next to the program name is the unique “Speed dial” number allocated to the program for rapid retrieval (see below).

## Programming Options

There are two options available for creating programs on the CYCLER:

1. Copy and edit one of the pre-written templates.
2. Create a new program from the defaults.

### Option 1: Creating programs using templates

There are two pre-written basic PCR templates: a two-step and a three-step template. These can be copied and modified as required. "In denat" stands for Initial Denaturing.

2 STEP		
Heated Lid	105°C	
Preheat Lid	on	
Pause	off	
- - - - -		
In denat	94.0°C	5m00
Hot Start	off	
- - - Stage 1 - - -		
Number cycles	30	
Step	94.0°C	0m30 max
Step	68.0°C	1m00 max
- - - - -		
Post run	72.0°C	5m00
Post hold	10.0 °C	

3 STEP		
Heated Lid	105°C	
Preheat Lid	on	
Pause	off	
- - - - -		
In denat	94.0°C	5m00
Hot Start	off	
- - - Stage 1 - - -		
Number cycles	30	
Step	94.0°C	0m30 max
Step	55.0°C	0m30 max
Step	72.0°C	0m30 max
- - - - -		
Post run	72.0°C	5m00
Post hold	10.0 °C	

Other templates available include Ice bucket (incubation at 4°C), Ligation (incubation at 15°C) and an RT-PCR template.

-*New*-		
2 STEP	[ 2 ]	<input type="checkbox"/>
3 STEP	3	<input type="checkbox"/>
[ ↑ ↓ , ← to select ]		

Use the "Down Arrow" key ↓ to select an existing 'Template' program from the menu, e.g. '2 Step Template' or '3 Step Template'.

Press the "Enter" key ↵ to accept it.

2 STEP	[ 2 ]
Run Method	
View Method	
[ ↑ ↓ , ← to select ]	

The Template is already defined and can be run as it is or easily copied, modified and saved under a new name.

```

      UNNAMED #
Heated Lid      105°C
Preheat Lid      on
[   ↑ ↓ , ← to edit   ]

```


A new name can now be given to the copied program (see below).

*To give a program a name*

```

■      UNNAMED #
Heated Lid      105°C
Preheat Lid      on
[   ↑ ↓ , ← to edit   ]


```

To type in a name for the program, press "Enter"  and a cursor will appear in the position shown.



Using the keypad, type in the name you wish to give the program.

The letters will scroll round with successive presses; for example the "2" key will give 'A' with one press; 'B' with two presses; 'C' with three presses and '2' with four presses. Five presses will again give 'A'.

When the name is complete, press  "Enter"

```

      DONNA
Heated Lid      105°C
Preheat Lid      on
[   ↑ ↓ , ← to edit   ]
[   ▽ to save         ]

```

There are now two options:


1. Save the program and run it as it is or
2. Edit the program then save and run it.

*To save the program*


```


      DONNA
Heated Lid      105°C
Preheat Lid      on
[   SAVE?       ]
[   ← YES or ▽ NO   ]

```

A program can only be run once it has been saved. Press the "End"  key.

The prompt line will flash between 'SAVE?' and 'YES or NO'.

Press "Enter"  for 'Yes' to save the changes you have made.



If you press "End"  for 'No' then all the changes you have made will be lost.

```

      DONNA      [7]
Run Program
View Program
[   ↑ ↓ , ← to select   ]

```

The program is now ready to run and will be allocated with a unique 'Speed dial' number which can be used to rapidly locate the program from the main start-up screen.

"Up Arrow" or "Down Arrow" keys   to scroll through the remaining list of options.



To edit a field or step in an existing program

**Methods** [ 0 ]  
 Information  
 Configuration  
 [ ↑ ↓ , ↵ to select ]

When the unit is switched on the display will show this screen.

If known, type in the 'Speed dial' number for the required program and press ↵ "Enter" This will take you directly to the required program.

DAVID 8 [D]  
 DONNA 7  
 [ ↑ ↓ , ↵ to select ]

If you do not know the 'Speed dial' number, press "Enter" ↵ to access the list of stored programs.

Either:

1. Scroll to the required program using the "Up Arrow" or "Down Arrow" keys or ↑ ↓
2. Type the first letter of the program name in the 'New program' field; this will then list all the program names beginning with that letter. Use the "Up Arrow" or "Down Arrow" keys ↑ ↓ to select the required program.

Press "Enter" to accept ↵

DONNA [ 7 ]  
 Run method  
 View method  
 Edit method  
 Copy method  
 Delete method  
 Set password  
 [ ↑ ↓ , ↵ to select ]

The full set of options for a program is shown here; these are shown on the LCD screen three lines at a time.

Use the "Down Arrow" and/or the "Up Arrow" ↑ ↓ keys to select the option you require.

Press the "Enter" key ↵ to select the option.

Edit a selected field or step

DONNA  
 Heated Lid 105°C  
 Preheat Lid on  
 [ ↑ ↓ , ↵ to edit ]  
 [ ▼ to save ]

Select 'Edit program' and a screen similar to this will appear.

Use the "Down arrow" and/or the "Up arrow" keys ↑ ↓ to select the field or step you want to edit.

You can now edit the active field or step by first pressing the "Enter" key ↵

```

DONNA
Heated Lid      ■ 105°C
Preheat lid      on
[ edit temperature ]

[ ← to disable ]

```

To edit a temperature, type in the required temperature. In the example here, the heated lid temperature is being edited.

Press “Enter” ↵ to accept.

Use the “Down Arrow” and/or the “Up Arrow” keys ↑ ↓ to select the next field or step you wish to edit.

```

DONNA
Heated Lid      105°C
Preheat lid      ■ on
[ ↑ ↓ to change ]

[ turn on/off ]

```

To edit a field with on/off settings, first select the field and press “Enter” ↵

Use the “Up Arrow” and “Down Arrow” ↑ ↓ keys to toggle between the settings.

Press “Enter” ↵ to accept.

```

Pause          off
-----
In denat  94.0°C  00m30
[ edit temperature ]

[ ← to disable ]

```

To edit a step with both a temperature and a time, first press “Enter” ↵ and type in the new temperature. Press “Enter” ↵ again. Now type in the new time. The default is minutes.

1. For whole minutes, type the number required then “Enter” ↵
2. To toggle between minutes and hours, use the “Decimal Point” key.
3. For values less than 1 minute, first press the “Decimal Point” key followed by the number of seconds required.
4. For minutes and seconds, type the number of minutes followed by the “Decimal Point” key then the number of seconds.

Once the values have been added, press “Enter” ↵ to accept.

```

Pause          off
-----
In denat  94.0°C  00m30
[ edit time ]

[ .for hrs/mins ]

```

Once you have finished editing and you are happy with the changes, press the “End” key ▽ followed by the “Enter” key ↵ to save.

```

-----
Post run  72.0°C  5m00
Post hold  10.0°C
[ ↑ ↓ , ↵ to edit ]

[ ▽ to save ]

```

If you press the “End” key ▽ a second time for ‘No’ then all the changes you have made will be lost.

To copy a selected program

```

      UNNAMED #
Heated Lid      105°C
Preheat Lid     on
[   ↑ ↓ , ← to edit   ]

[   ▽ to save         ]

```

To delete a selected program

To delete a selected program

```

      DONNA

delete program
SURE?

[   ← Yes or ▽ No     ]

```

To set a password for a selected program

```


      DONNA

New passwd  █
[   ← to accept       ]


[   ▽ to exit         ]

```


If you select 'Copy program' a screen similar to this will appear and the software will automatically give the program the next number until it is named.

Press "Enter"  and type in a new name. You will then be able to edit the program as previously described.

If you select 'Delete program' a screen similar to this will appear.

Press "Enter"  to delete the program or "End"  to return to the program screen.

If you want to password protect the program so that other users cannot edit it, select 'Set password'. A screen similar to this will appear.

Type in the password you wish to give it and press "Enter" . The program list will then display the key symbol next to the name to show that it is locked.

Other users will still be able to copy/view your program but not edit it.

## Option 2: Creating a new program

A second option is to create a completely new program without using templates.

```

      UNNAMED #
Heated lid      105°C
Preheat lid     off
Pause          off
- - - - -
In denat      94.0°C 5m00
Hot Start      off
- - - - -
Post run      72.0°C 5m00
Post hold      10.0 °C

```

A number of typical values are pre-entered including:

Heated lid temperature

Preheat lid

Initial denaturation

Hot start

Final extension

Final hold

These values can be edited and the additional stages and steps required for cycling can be added by following the instructions detailed on the following pages.

## Example programs

A couple of example thermal cycling programs are shown below.

### Typical 3-Step PCR

```
Heated Lid      105
Preheat Lid     on
Pause           off
-----
In denat   94°C      5min
Hot Start                      off
----- Stage 1 -----
Number Cycles                      30
Step       94°C      0m30      max
Step       55°C      0m30      max
Step       72°C      0m30      max
-----
Post run           72°C      5m
Post hold          4°C      10m
```

### 3-Step PCR with multiple 'Stages'

```
Heated Lid      105
Preheat Lid     on
Pause           off
-----
In denat   94°C      5min
Hot Start                      off
----- Stage 1 -----
Number Cycles                      10
step       94°C      0m30      max
step       60°C      0m30      max
step       72°C      0m30      max
----- Stage 2 -----
Number Cycles                      20
step       94°C      0m30      max
step       55°C      0m30      max
step       72°C      0m30      max
-----
Post run           72°C      5m
Post hold          4°C      10m
```

## Programming

```
- *New* - [ ]
2 STEP   2
3 STEP   3
[↑ ↓ , ← to select ]
```

Create a new program by selecting  
-\* New \*- from the menu and press "Enter"



```
UNNAMED #
Heated Lid      105°C
Preheat Lid     on
[ ↑ ↓ , ← to edit ]
[ ▼ to save ]
```

Give the program a name and edit any of  
the default fields as described above.

To add a stage and then edit it

```
In denat   94.0°C      5m00
Hot Start                      off
-----
[ ← to insert stage ]
```

Use the "Down Arrow" ↓ key to scroll to  
the dotted line below 'Hot Start'; it will  
start to flash.

Press "Enter" ← to insert a stage

```

----- Stage 1 -----
Number cycles          █
Step  94.0°C 00m30  max
[    ← edit no.cycles  ]

```

A new stage will be inserted which consists of a 'Number cycles' field and one step.

Edit the number of cycles by typing the number required then press "Enter"↵

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[    '.' to insert step  ]

[    ↑↓, ← to edit      ]
[    ← to delete step   ]

```

Use the "Down Arrow"↓ key to move to the step.

Press "Enter"↵ to edit the step as described above.

Press "Enter"↵ again to accept.

---

#### *To insert additional steps*

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[    '.' to insert step  ]

[    ↑↓, ← to edit      ]

```

To insert a new step before the first step, use the "Up Arrow" or "Down Arrow"↑ ↓ keys to select 'Number cycles'.

Press the "Decimal Point" key to insert a new step.

To insert a new step after an existing step, use the "Up Arrow" or "Down Arrow"↑ ↓ keys to select the step.

Press the "Decimal Point" key to insert a new step.

---

#### *To delete a step*

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[    '.' to insert step  ]

[    ↑↓, ← to edit      ]
[    ← to delete step   ]

```

Use the "Up Arrow" or "Down Arrow"↑ ↓ keys to select the step you wish to delete.

Press "Delete"← to delete the step.

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[    DEL STEP ?        ]

[    ← Yes or ▽ No     ]

```

The prompt line will flash between 'DEL STEP?' and 'YES or NO'.

Press "Enter"↵ to delete the step.  
or

Press "End"▽ to return to editing the step.

*To add a new stage before an existing one*

---

```
----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[  ← to insert stage  ]

[  ← to delete stage  ]
```

Use the "Up Arrow" or "Down Arrow" keys to highlight the stage number. ↑ ↓

Press "Enter"↵

A new stage will be added before the existing stage. The other stage numbers will be adjusted as appropriate. The new stage can now be edited as described above.

*To add a new stage at the end*

---

```
- - - - -
Post run 72.0°C  5m00
Post hold      10.0°C
[  ↵ to insert stage  ]
```

Use the "Up Arrow" or "Down Arrow"↑ ↓ keys to highlight the dotted line below the last stage.

Press "Enter"↵

A new stage will be added with the next consecutive number. The new stage can now be edited as described above.

*To delete a stage*

---

```
----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[  ← to insert stage  ]

[  ← to delete stage  ]
```

Use the "Up Arrow" or "Down Arrow"↑ ↓ keys to highlight the stage you wish to delete.

Press "Delete"←

```
----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[      DEL STAGE?      ]

[  ↵ Yes or ▽ No  ]
```

The prompt line will flash between 'DEL STAGE?' and 'YES' or 'NO'.

Press "Enter"↵ to delete the stage.

*To insert a 'pause' into a program*

---

```
Step  94.0°C  00m30  max
----- Stage 2 -----
Number cycles          █
Step  94.0°C  00m30  max

[  '.' to insert step  ]
[  ↑ ↓ , ← to edit   ]
```

To insert a pause after an existing stage, follow the steps for inserting a stage as above.

Highlight the 'Number cycles' line then press "Pause" **||**

```
Step  94.0°C  00m30  max
----- Stage 2 -----
Number cycles          0
Step  94.0°C  00m30  max

[  CHANGE TO PAUSE    ]
[  ← Yes or ▽ No     ]
```

The prompt line will flash between 'CHANGE TO PAUSE?' and 'YES or NO'.

Press "Enter" **↵** to insert a pause.


```
----- Stage 2 -----
Pause
-----
[  ← to insert step  ]

[  ← to delete step  ]
```

The stage number will then be highlighted with 'Pause' as the only line.


## To program increment or decrement temperature or time.

Under normal circumstances, the Hold Temperature of all steps is constant. However, it is possible to automatically increment or decrement the temperature of a specified step of a programme.

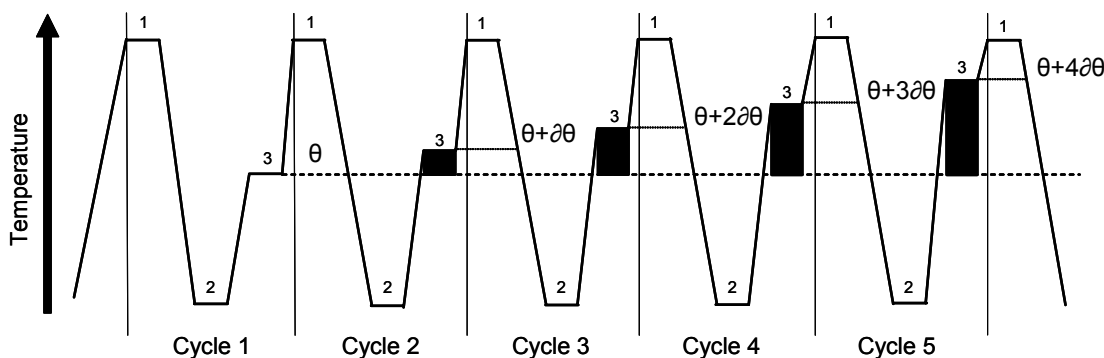

*Temperature profile of incremented Hold Temperature at step 3 of each cycle*

The Hold Temperature of the incremented/decremented step is the defined Hold Temperature plus the summation of the increments/decrements. The first cycle is never incremented/decremented, only subsequent cycles. If you select a decrement, the Hold Temperature is prevented from falling below 4°C. If you select an increment, the Hold Temperature is prevented from rising above 99°C. Similarly, under normal circumstances, the Hold Time of all steps is constant. However, it is possible to automatically increment or decrement the duration of a specified step of a program.

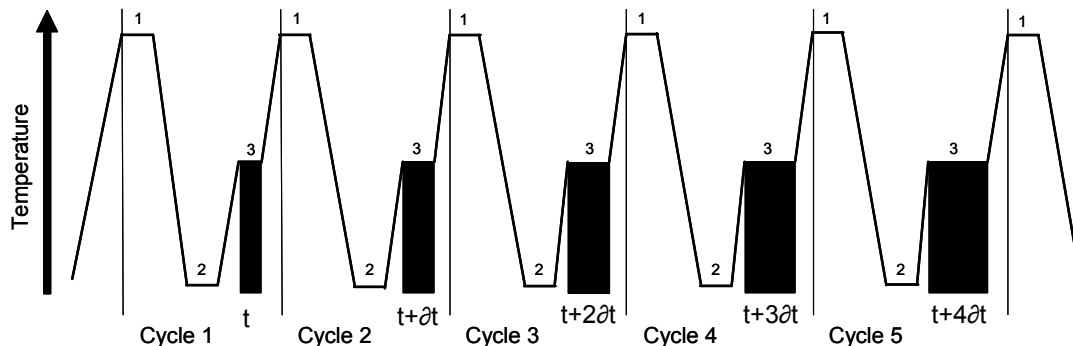

*Temperature profile of incremented Hold Time at step 3 of each cycle*

The Hold Time of the incremented/decremented step is the defined Hold Time plus the summation of the increments/decrements. The first cycle is never incremented/decremented, only subsequent cycles. If you select a decrement, the Hold Time is prevented from falling below one second.

Number cycles	30
Step 94.0°C	00m30 max
Step 58.0°C	00m30 max
[ , to edit ]	

[ ← to delete step ]  
[ '.' to insert step ]

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to move to the step where you want an increment or decrement temperature or time.

Press "Enter" ↵ to go into edit mode.

Press "Pause" ||



```

Number cycles          30
Step  94.0°C   00m30   max
Step  58.0°C   00m30   max
[    CHANGE STEP ?    ]

[    ↵ Yes or ▽ No    ]

```

The prompt line will change to flash between 'CHANGE STEP?' and 'YES' or 'NO'.

Press "Enter"↵

The cursor will again move into the edit mode.

Now press the "Up Arrow" key ↑

```

Number cycles          30
Step  94.0°C   00m30   max
fst   55.0°C   00m30   max
[    ↑ ↓, ↵ to edit    ]

[    ▽ to save         ]

```

The step will change to 'fst', meaning "first", and the software will insert another line after it entitled 'lst', meaning "last".

Press "Enter"↵ twice to go into edit mode.

```

Number cycles          30
Step  94.0°C   00m30   max
fst   65.0°C   00m30   max
[    edit temperature  ]

```

Use the "Alpha/numeric" keys to change the values of the 'fst' line to the first temperatures and/or times you require.

Press "Enter"↵ after each field you change.

You will exit 'Edit Mode' when you reach the end of the line.

```

Step  94.0°C   00m30   max
fst   65.0°C   00m30   max
lst   55.0°C   00m30
[    ↑ ↓, ↵ to edit    ]

[    ← to delete step  ]
[    \.' to insert step ]

```

Use the "Down Arrow" key ↓ to move onto the 'lst' line.

Press "Enter"↵ to edit the step.

```

Step  94.0°C   00m30   max
fst   65.0°C   00m30   max
lst   55.0°C   1m30
[    edit time         ]

[    .For hrs/mins     ]

```

Use the "Alpha/numeric" keys to change the values of the '1st' line to the final temperatures and/or times you require.

Press "Enter"↵ for each field you change.

You will exit 'Edit Mode' when you reach the end of the line.

When you run the program it will automatically increment or decrement in equal steps from the first time and/or temperature to the last over the number of cycles in the stage. For example: if the

first temperature is 65°C and the last temperature is 55°C, over 30 cycles the temperature will decrease 0.33°C each cycle. Similarly, in the example shown, the time would increase from 30 seconds to 1 minute 30 seconds in 2 second increments.

---

*To save the program*

---

```
----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[                SAVE?                ]

[  ← Yes or ▽ No  ]
```

A program can only be run once it has been saved. Press the ▽ “End” key.

The prompt line will flash between 'SAVE?' and 'YES or NO'.

Press “Enter” ↵ for ‘Yes’ to save the changes you have made.

If you press “End” ▽ for ‘No’ then all the changes you have made will be lost.

```
GEORGE [10]
Run method
View method
[↑ ↓ , ↵ to select ]
```

The program is now ready to run and will be allocated with a unique “Speed dial” number which can be used to rapidly locate the program from the main start-up screen.

Use the “Up Arrow” and “Down Arrow” ↑ ↓ keys to scroll through the list of options.

---

*Stopping what you are doing*

---

At any time during programming the CYCLER, when you are satisfied with the program, you can finish and accept what you have done by first pressing the “End” key. ▽

Then follow the instructions on screen. If you are not sure, wait for the instructions in the prompt line to scroll back to the first line that you see.

## To run a program

```
Methods [0]
Information
Configuration
[↑ ↓ , ← to select ]
```

When the unit is switched on the display will show this screen.

```
Methods [7]
Information
Configuration
[↑ ↓ , ← to select ]
```

If known, type in the 'Speed dial' number for the required program and press ← "Enter" This will take you directly to the required program.

If you do not know the 'Speed dial' number, press "Enter" ← to access the list of stored programs.

```
* * New Methods * * [ ]
2 STEP 2
3 STEP 3
[↑ ↓ , ← to select ]
```

Next either:

1. Scroll down to the required program using the "Up Arrow" and "Down Arrow" ↑ ↓ keys.
2. Or type the first letter of the program name in the "New program" field; this will then list all the program names beginning with that letter. Use the "Up Arrow" and "Down Arrow" keys ↑ ↓ to select the required program.

```
DAVID 8 [D]
DONNA 7
[ ↑ ↓ , ← to select ]
```

Press "Enter" to accept ↓  
For example 'Donna'.

```
DONNA [7]
Run method
View method
[↑ ↓ , ← to select ]
```

To run the program, select 'Run program' and press "Enter" ↓

```
DONNA
[ RUN METHOD ? ]
[ ← Yes or ▽ No ]
```

A screen similar to this one will appear.

Press "Enter" ↓ to start the run.

```

Cycle 4/30      56m08
Sample 58.0°C   5.7□ /S
Hold @ 58.0°C   00m17
[  ▽ to STOP   ]
[  || to PAUSE  ]
[  ↵ for info   ]

```

During the run the LCD screen will report:

1. The cycle number
2. The time remaining (counting down in seconds)
3. The sample temperature (whether holding or ramping)
4. Maximal ramp rate
5. The current step of the run
6. A count down of the hold time.

An example is shown here.

---

### Manual pause

```

Cycle 24/30      19m26
Sample 72.0°C    2.7□ /S
Hold @ 72.0°C    00m10
[  PAUSE ?      ]
[  ↵ YES or ▽ NO ]

```

To pause a program while it is running, press "Pause" **||**

The run will continue but the prompt line will alternate 'PAUSE?' and 'YES or NO'

To confirm that you want the program to pause press "Enter" **↵**

If you have pressed the "Pause" key by mistake and you want the program to continue, press "End" **▽**

The unit will emit a beeping sound while paused.

To resume the run, press "Pause" **||** followed by "Enter" **↵** to continue.

```

Cycle 24/30      19m26
Sample 72.0°C    5.7□ /S
Paused
[  ▽ to STOP     ]
[  || to CONTINUE ]
[  ↵ for info     ]

```

---

### Manual Stop

```

Cycle 28/30      10m39
Sample 94.0°C    5.7□ /S
Hold @ 94.0°C    00m10
[  STOP PROGRAM? ]
[  ↵ YES or ▽ NO ]

```

To stop a program while it is running, press "End" **▽**

The run will continue but the prompt line will alternate 'STOP PROGRAM?' and 'YES or NO'

To confirm that you want to stop the program press "Enter" **↵**

If you have pressed the "End" **▽** key by mistake and you want the program to continue, press "End" **▽** again.

When a program is stopped the unit will emit a beeping sound and the original program screen will appear.

```
          DONNA
Heated Lid           on
Status               ramping
[      ↵ to exit    ]
```

Pressing “Enter” ↵ while a program is running will give information about the current run including which program is actually running and the unit status e.g. holding or ramping.

An example is shown here.

## 7. Information

```
Methods              [0]
Information
Configuration
[↑ ↓ , ↵ to select ]
```

The Information field on the front screen contains all the unit-specific information such as serial number and software version.

Press “Enter” ↵ to access the unit information.

```
Block Type:      48x0.2ml
Block id         123456789
Unit s/n         C3C123456
S/w version      6.zz
H/w version      2.ZZ
Comms address    1
Unit cycles      624
Block cycles     69
```

The complete list is displayed here.

## 8. Configuration

**Warning!** The Configuration software should only be accessed by designated personnel, as some parameters can be changed or may be lost when using it.

```
Methods              [0]
Information
Configuration
[↑ ↓ , ↵ to select ]
```

The Configuration field on the front screen allows a system administrator access to set up all the default settings for the programs, time and date settings and to change and remove passwords.

```

CONFIGURATION
Password 1234
[  ← to accept  ]

```

```

[  ▽ to exit  ]

```

```

Settings
Passwords
Format memory
Self-calibrate

```

Access to the 'Configuration' set up is password protected.  
 Default admin passwords is 1234  
 To enter configuration, use the alpha keys to type in '1234'

Press "Enter" ↵

There are a number of options which can be changed; these are listed here.

Use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to select then press "Enter" ↵ to access any of the options. Each option is discussed in detail below.

### Settings

---

```

Settings
Passwords
Format memory
Self-calibrate

```

Highlight 'Settings' then press "Enter" ↵

```

Settings
Passwords
Format memory
Self-Calibrate

```

Highlight 'Settings' then press "Enter" ↵

```

Set date      xx/xx/xx
Set time      xx:xx
Auto restart  on
Beeper        on
Coms address  1
Coms speed    19200
[↑ ↓ , ← to select ]

```

The complete list of settings is given here.

Auto restart "on" will allow the unit to restart from where it left off should there be a power failure and the power returns.

The "Comms address" refers to the location of a unit in a daisy chain and must be unique if there are several units linked to each other.

The optimum "Coms speed" is set at 19200. When using a control software, ensure all cyclers are set to the same value.

Edit any line as described for 'Program defaults' above. Press "Enter" ↵ to accept the new value.

When you have finished, press the “Enter” key to go back to the ‘Configuration’ menu ↵

## Passwords

```
Settings
Passwords
Format memory
Self-calibrate
```

The ‘Passwords’ function allows the administrator to change the ‘Admin password’ (which allows access to Configuration) and also to delete any passwords associated with programs.

Highlight ‘Passwords’ then press “Enter” ↵

The screen will appear as opposite.

We suggest that you change the ‘Admin password’ at this stage in order to protect the configuration settings.

```
* * Admin passwd * *
DONNA
GEORGE
[ ↑ ↓ , ← to select ]
```

```
ADMIN PASSWD
Password █
Confirm
[ ← to accept ]
```

Use the “Up Arrow” ↑ key to highlight ‘Admin passwd’

Press “Enter” ↵ and type in the new password. Repeat to confirm and accept.

The screen will go back to the ‘Configuration’ menu.

```
* * Admin passwd * *
DONNA
GEORGE
[ DEL PASSWORD? ]
[ ← Yes or ▽ No ]
```

To delete a password associated with a program, use the “Up Arrow” and “Down Arrow” ↑ ↓ keys to highlight the program.

Press “Enter” ↵

Press “Enter” ↵ again to delete the password.

## Controller reset

```
Settings
Passwords
Format memory
Self-calibrate
```

The ‘Format memory’ will delete all saved programs from the unit.

Highlight ‘Format memory’ then press ↵ “Enter”

```
Format Memory
This will delete
all methods
[ SURE ? ]
[ ← YES ▽ NO ]
```

To delete all the programs, press “Enter” ↵  
Set ADMIN PASSWORD to 1234.

## 9. Notes on transferring protocols from other thermal cyclers

The CYCLER TCY heats at over 3.6°C/s and provides the highest levels of temperature uniformity. Many slower machines have historically begun the countdown on hold times before the set temperature is reached in order to avoid any overshooting in sample temperature. The slower the heating rate the more significant the loss in hold time, i.e. if a 30 second hold time is set at 55°C then perhaps 5 to 10s will be lost in between the start of the time countdown and actually reaching the set temperature. The overall effect of this method of control is that your sample, if cooling from a higher temperature, is subjected to an environment with an elevated temperature compared to the target temperature for a significant part of the hold time.

The CYCLER TCY provides the user with 'what-you-program-is what-you-get'. Due to the rapid and responsive nature of the Peltier heating system, the time between the timer countdown initiating and actually reaching the set temperature has been reduced to as little as 1s. This ensures that the sample is only subjected to the exact temperature required.

Taking the above into consideration it must therefore be noted that care must be taken when re-optimizing protocols from a slower machine.

### Trouble Shooting

1. The optimum annealing temperature may in fact be higher than the set temperature used in the slower machine. Try increasing it by 1 or 2°C.
2. The hold time can be reduced once the optimum temperature has been found. Try reducing it by 5 to 10s.



## 10. User maintenance


### **Cleaning your CYCLER**

The heating/cooling block, including wells and flat surfaces, should be cleaned regularly to ensure optimum heat transfer to the samples. Always clean the block if there has been a spillage. Use a cloth or cotton buds dipped in a fresh, 50:50 water/isopropanol solution and make sure that no deposits are left in the wells.

In the case of radioactive spillages the heater block can be removed from the unit for more detailed cleaning. Omega Engineering, Inc. recommend that you use a proprietary cleaning agent and follow the manufacturer's instructions. The heating/cooling block is made of aluminum, therefore an agent such as Neutracon (from Decon Labs, Inc.), suitable for nonferrous metals should be used. However, remember that other parts of the unit are made of ferrous materials and may be damaged by spillage onto them.

### Heater Block removal

Before cleaning your unit, disconnect from the power supply and allow it to cool to below 50°C. Remove the four screws on the underside of the unit and drop the block/fan assembly down. Unplug the assembly from the wiring and remove the assembly from the unit. Refitting the block/fan assembly is the reverse procedure; be careful of the wires as you refit the block.

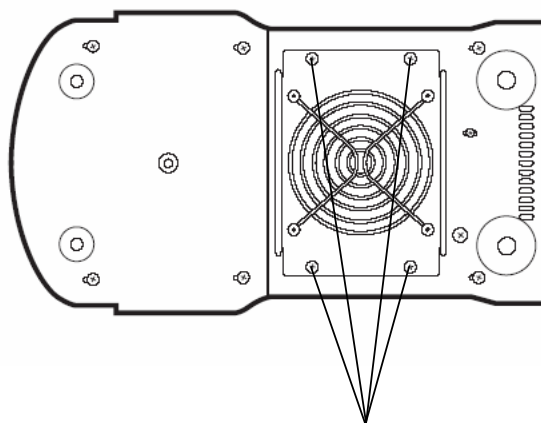

***Underside of the CYCLER TCY***

Four screws which hold  
the block/fan assembly in place

The outer case of the CYCLER TCY can be cleaned with a cloth dipped in water or ethanol (methanol or formaldehyde can also be used). No part of the case or cover should be immersed in the solvents. Do not use aggressive solvents such as acetone, or abrasive cleaners.

## 11. Some frequently asked questions

- Q1** What is the 'Pause' function at the start of the program used for?
- A1** Some users prefer to preheat the heated lid before placing the samples into the unit. The pause feature is used to stop the unit after the 4 minute heated lid preheat step. It will also sound an audible alarm indicating that the machine is ready for the sample tubes or plate to be added. Pressing the "Pause" key followed by the "Enter" key will commence the remaining program.
- Q2** Why do you have to press some keys twice to make the unit do what you want it to do?
- A2** With many of the program features there is a fail-safe function which asks if you are sure you want to perform a particular action. You are required to press one key to select a function followed by a second to accept it.
- Q3** What is 'Hot Start'?
- A3** A "hot start" is used for reducing non-specific product formation during the first cycle of amplification. The 'Hot Start' programming step is used to pause the instrument at a specific temperature, typically around 70°C, after the initial template denaturation. The reason is to allow the manual addition of unmodified *Taq* DNA polymerase which may lose activity if added during the initial 5min denaturation. Heat-activated *Taq* or Hot Start enzymes do not require this step.
- Q4** What is the incremental function? (See page 24).
- A4** Incremental timing and temperature are used to increase or decrease either the time or temperature incrementally over the number of cycles in a stage. Incrementation of extension time is used with 'Long PCR' which is when large template fragments are to be amplified (e.g. 27 kb lambda DNA, 40kb genomic DNA).  
Decremental temperature is used for protocols such as 'Touchdown PCR' where one starts with a high annealing temperature in the first cycle and gradually decreases the temperature over the number of cycles in the stage. This ensures that only the specific product is amplified.
- Q5** The heat-up ramp-rate in the specification is 3.6°C/second yet I can only manually set it to 2.0°C/second. Why is this?
- A5** If you set the ramp-rate to max the block will heat at 3.6°C/second the unit sets its own control parameters. If you program a ramp-rate the control parameters reduce the ramp-rate. This is a feature of the software and not the actual ramp-rate.

## **12. Additional information**

Brief fault finding notes and a list of replacement parts are given in this section.

Note that this equipment should only be dismantled by properly trained personnel. Removing the outer cover exposes potentially lethal power voltages.

There are no user serviceable parts within this equipment.

### **Fault Finding**

Should you encounter problems with your TCY contact Omega Flow Engineering for assistance.

### **Fuses**

If neither the power light nor display on the front panel is lit, one of the two fuses may have blown. Check that there is no external cause, such as a faulty plug or lead. Check both fuses and replace the faulty fuse with a new one of the correct value (fuse values are given on the label next to the power inlet). Note that fuses should only be replaced by a qualified electrician.

The holder for the two fuses is built into the power input socket. First remove the power cable and then gently prise the fuse drawer open with a flat-bladed screwdriver or similar tool.

Each fuse can be removed by using the screwdriver as a lever.

Exchange the faulty fuse in the fuse holder for a working fuse of the correct value. Finally, replace the fuse drawer in the fuse compartment and push the drawer shut.

Fuses which blow repeatedly indicate a serious fault and you should return the unit to your supplier for repair.

### **The heated lid over-temperature cut-out**

The heated lid is fitted with an independent circuit to protect it from overheating. In the unlikely event of an over-temperature problem with the lid, the unit is fitted with a thermal fuse which will remove power to the heater plate should the maximum temperature be exceeded.

### **Insulation testing**

This equipment is fitted with RFI suppression circuitry. Any check of the electrical insulation by means of high voltage dielectric testing (for example as in BS EN 61010-1) must be carried out using only a TCY voltage.

This unit contains semiconductor components which may be damaged by electric field effects.

### **Interchangeable blocks**

The block can be removed and replaced by another of the same sort or of a different sort. The software will register the change and set up the unit for the new block.

Remove the four screws on the underside of the unit and drop the block/fan assembly down. Unplug the assembly from the wiring and remove the assembly from the unit. Refitting the block/fan assembly is the reverse procedure; be careful of the wires as you refit the block.

## WARRANTY/DISCLAIMER

OMEGA ENGINEERING, INC. warrants this unit to be free of defects in materials and workmanship for a period of **36 months** from date of purchase. OMEGA's WARRANTY adds an additional one (1) month grace period to normal **three (3) year product warranty** to cover handling and shipping time. This ensures that OMEGA's customers receive maximum coverage on each product.

If the unit malfunctions, it must be returned to the factory for evaluation. OMEGA's Customer Service Department will issue an Authorized Return (AR) number immediately upon phone or written request. Upon examination by OMEGA, if the unit is found to be defective, it will be repaired or replaced at no charge. OMEGA's WARRANTY does not apply to defects resulting from any action of the purchaser, including but not limited to mishandling, improper interfacing, operation outside of design limits, improper repair, or unauthorized modification. This WARRANTY is VOID if the unit shows evidence of having been tampered with or shows evidence of having been damaged as a result of excessive corrosion; or current, heat, moisture or vibration; improper specification; misapplication; misuse or other operating conditions outside of OMEGA's control. Components in which wear is not warranted, include but are not limited to contact points, fuses, and triacs.

**OMEGA is pleased to offer suggestions on the use of its various products. However, OMEGA neither assumes responsibility for any omissions or errors nor assumes liability for any damages that result from the use of its products in accordance with information provided by OMEGA, either verbal or written. OMEGA warrants only that the parts manufactured by the company will be as specified and free of defects. OMEGA MAKES NO OTHER WARRANTIES OR REPRESENTATIONS OF ANY KIND WHATSOEVER, EXPRESSED OR IMPLIED, EXCEPT THAT OF TITLE, AND ALL IMPLIED WARRANTIES INCLUDING ANY WARRANTY OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE ARE HEREBY DISCLAIMED. LIMITATION OF LIABILITY: The remedies of purchaser set forth herein are exclusive, and the total liability of OMEGA with respect to this order, whether based on contract, warranty, negligence, indemnification, strict liability or otherwise, shall not exceed the purchase price of the component upon which liability is based. In no event shall OMEGA be liable for consequential, incidental or special damages.**

CONDITIONS: Equipment sold by OMEGA is not intended to be used, nor shall it be used: (1) as a "Basic Component" under 10 CFR 21 (NRC), used in or with any nuclear installation or activity; or (2) in medical applications or used on humans. Should any Product(s) be used in or with any nuclear installation or activity, medical application, used on humans, or misused in any way, OMEGA assumes no responsibility as set forth in our basic WARRANTY/DISCLAIMER language, and, additionally, purchaser will indemnify OMEGA and hold OMEGA harmless from any liability or damage whatsoever arising out of the use of the Product(s) in such a manner.

## RETURN REQUESTS/INQUIRIES

Direct all warranty and repair requests/inquiries to the OMEGA Customer Service Department. BEFORE RETURNING ANY PRODUCT(S) TO OMEGA, PURCHASER MUST OBTAIN AN AUTHORIZED RETURN (AR) NUMBER FROM OMEGA'S CUSTOMER SERVICE DEPARTMENT (IN ORDER TO AVOID PROCESSING DELAYS). The assigned AR number should then be marked on the outside of the return package and on any correspondence.

The purchaser is responsible for shipping charges, freight, insurance and proper packaging to prevent breakage in transit.

FOR **WARRANTY** RETURNS, please have the following information available BEFORE contacting OMEGA:

1. Purchase Order number under which the product was PURCHASED,
2. Model and serial number of the product under warranty, and
3. Repair instructions and/or specific problems relative to the product.

FOR **NON-WARRANTY** REPAIRS, consult OMEGA for current repair charges. Have the following information available BEFORE contacting OMEGA:

1. Purchase Order number to cover the COST of the repair,
2. Model and serial number of the product, and
3. Repair instructions and/or specific problems relative to the product.

OMEGA's policy is to make running changes, not model changes, whenever an improvement is possible. This affords our customers the latest in technology and engineering.

OMEGA is a registered trademark of OMEGA ENGINEERING, INC.

© Copyright 2009 OMEGA ENGINEERING, INC. All rights reserved. This document may not be copied, photocopied, reproduced, translated, or reduced to any electronic medium or machine-readable form, in whole or in part, without the prior written consent of OMEGA ENGINEERING, INC.

# Where Do I Find Everything I Need for Process Measurement and Control? **OMEGA...Of Course!**

*Shop online at [omega.com](http://omega.com)<sup>SM</sup>*

## **TEMPERATURE**

- ☑ Thermocouple, RTD & Thermistor Probes, Connectors, Panels & Assemblies
- ☑ Wire: Thermocouple, RTD & Thermistor
- ☑ Calibrators & Ice Point References
- ☑ Recorders, Controllers & Process Monitors
- ☑ Infrared Pyrometers

## **PRESSURE, STRAIN AND FORCE**

- ☑ Transducers & Strain Gages
- ☑ Load Cells & Pressure Gages
- ☑ Displacement Transducers
- ☑ Instrumentation & Accessories

## **FLOW/LEVEL**

- ☑ Rotameters, Gas Mass Flowmeters & Flow Computers
- ☑ Air Velocity Indicators
- ☑ Turbine/Paddlewheel Systems
- ☑ Totalizers & Batch Controllers

## **pH/CONDUCTIVITY**

- ☑ pH Electrodes, Testers & Accessories
- ☑ Benchtop/Laboratory Meters
- ☑ Controllers, Calibrators, Simulators & Pumps
- ☑ Industrial pH & Conductivity Equipment

## **DATA ACQUISITION**

- ☑ Data Acquisition & Engineering Software
- ☑ Communications-Based Acquisition Systems
- ☑ Plug-in Cards for Apple, IBM & Compatibles
- ☑ Datalogging Systems
- ☑ Recorders, Printers & Plotters

## **HEATERS**

- ☑ Heating Cable
- ☑ Cartridge & Strip Heaters
- ☑ Immersion & Band Heaters
- ☑ Flexible Heaters
- ☑ Laboratory Heaters

## **ENVIRONMENTAL MONITORING AND CONTROL**

- ☑ Metering & Control Instrumentation
- ☑ Refractometers
- ☑ Pumps & Tubing
- ☑ Air, Soil & Water Monitors
- ☑ Industrial Water & Wastewater Treatment
- ☑ pH, Conductivity & Dissolved Oxygen Instruments